

Membrane filtration – mastering microbiology and improving asset utilisation with BactoSense



Abstract

Filtration systems, such as ultrafiltration and reverse osmosis, are critical in producing clean water by removing particulate matter and bacteria. However, their performance can be compromised by operational conditions, biofouling and membrane breaches, requiring continuous monitoring to ensure consistent water quality at the point of use. Traditional methods, like heterotrophic plate count (HPC), are time consuming and offer delayed insights. In contrast, the BactoSense solution provides real-time bacterial monitoring with precise and actionable results in just 20 minutes. In this application note, we outline key application areas and accomplishments reported by BactoSense users in drinking water production, industrial water systems, water reuse and desalination.

Keywords

End-use: drinking water, water reuse, desalination, industrial water
Process: microfiltration, ultrafiltration, nanofiltration, reverse osmosis

Filtration for microbiological purification

Filtration systems are commonly used to remove particulate matter from water. In terms of microbiological purification, membrane filtration methods like ultrafiltration and reverse osmosis effectively eliminate bacteria, while media filters such as activated carbon filters help by removing nutrients that could otherwise promote microbial growth. As a result, filtration systems play a crucial role in producing clean and safe water.

However, filtration systems are not foolproof. Their effectiveness depends on a complex interaction of factors, including operational conditions, the properties of the feed water, defects, breaches, biofouling and maintenance. Continuous monitoring is essential to consistently meet water quality standards at the point of use. Traditional growth-based methods, such as heterotrophic plate count (HPC), must be conducted in a lab, are tedious and take several days. In contrast, with BactoSense, bacteria are directly counted and precise results are available in just 20 minutes.

BactoSense to increase asset utilisation

BactoSense has been developed to make microbiological data actionable. Imagine having a tool in your hands to continuously verify the bacterial removal (e.g. log reduction values) of a membrane system, immediately detect filter breaches or declining filtrate quality, swiftly qualify individual assets, and gain a complete understanding of the impact of operational changes on the microbiological quality of the filtrate.

The benefits of BactoSense have been demonstrated across various applications, from monitoring membrane integrity to optimising operational parameters and maintenance practices. Key application areas and accomplishments reported by BactoSense users include:

Filtrate monitoring

Online bacterial monitoring with measurement intervals as short as 30 minutes allows operators to be alerted immediately if filtrate quality is abnormal.

- Verify the integrity of membrane filtration systems
- Implement an early warning system to detect membrane breaches and gradual membrane performance loss, e.g. due to biofouling or scaling.
- Ensure the proper functioning of filtration barriers under all operational conditions, e.g. during maintenance activities or other upstream operations.

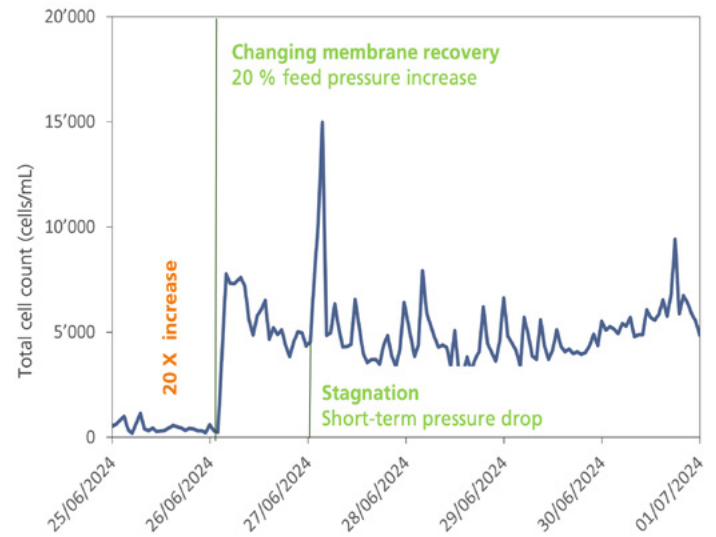


Figure 1: Measuring reverse osmosis permeate quality during process optimisation. To increase membrane recovery, the feed pressure was increased by 20%, resulting in an immediate 20-fold increase in the total cell count. Importantly, this phenomenon was not transient. The cell concentrations remained at this high level, indicating that operational parameters can strongly influence the microbiological quality of the permeate. In addition, short-term stagnation was shown to lead to bacterial regrowth and lowered permeate quality.

Process optimisation

First, understand your process comprehensively, and then improve its performance and resilience.

- Optimise timings for filter backwashes and duration of subsequent discharge.
- Investigate the effect of operational changes (e.g. flow rate or recovery) on the microbiological filtrate quality.
- Understand the kinetics of microbial regrowth during stagnation.
- Quantify the filtration efficiency under all relevant operational conditions by measuring feed and filtrate.

Troubleshooting

There is no need to wait days for lab results. Getting results in only 20 minutes speeds up your troubleshooting and helps reduce downtime, whether offline, atline, or online.

- After detecting an anomaly in a combined filtrate, directly investigate individual membrane modules.
- Swiftly identify poorly performing or defective membranes to mitigate the risk of pathogen passage.
- Locate the potential source of contamination, such as membrane failure, regrowth in the feed storage tank, or biofouling of a prefilter.

Asset commissioning, qualification and life cycle management

Things can go wrong during the commissioning of new assets. Implement a monitoring programme to make sure you are in control of microbiology.

- Qualify your filtration system: for membrane filtration systems, determine log reduction values by measuring feed and filtrate concentrations online or at-line. Ensure performance requirements are met.
- Continuously monitor the filtrate of newly installed membranes to detect “burn-in” failures swiftly.

Feed and filtrate

Better understand the properties of feed and filtrate to optimise filter operation and ensure filter performance requirements are met under all conditions.

- Assess the microbiological fluctuations in the feed water to improve filter operation (e.g. seasonal or operational variation).
- Locate microbial contaminations in feed tanks or bio-fouling in prefilters.
- Understand and manage microbial regrowth downstream until reaching the point of use.
- Assess the biostability of the water in various steps of the treatment process.

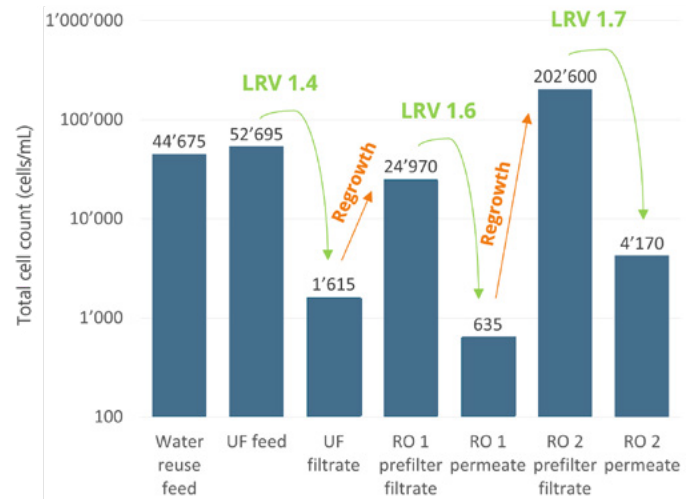


Figure 2: Microbiological characterisation of a water reuse treatment train, consisting of ultrafiltration followed by two-stage reverse osmosis. Membrane treatment steps significantly reduced total cell counts, but subsequent microbiological regrowth in permeate tanks and reverse osmosis prefilters drastically increased bacterial levels again, lowering the overall bacterial log reduction across the treatment train. LRV, log reduction value; RO, reverse osmosis; UF, ultrafiltration.

The examples above demonstrate that BactoSense is a game changer for filtration system monitoring and management. Unlike traditional methods, it empowers operators to detect membrane issues early, optimise performance and swiftly address potential risks. By providing continuous, actionable insights, BactoSense revolutionises how water quality is managed in processes such as drinking water production, desalination and water reuse.

Safe Water. Anytime. Anywhere.

bNovate Technologies SA

Ch. Dent d'Oche 1A · CH-1024 Ecublens
Tel. +41 (0)21 552 14 21
info@bnovate.com · www.bnovate.com